


Saliva Collection Instructions

Do NOT eat, drink, smoke, or chew gum for 30 minutes before giving your saliva sample.



Fill the tube with saliva to the black wavy line.

Fill the tube until your saliva (not including bubbles) is at or just above the wavy line. **DO NOT OVERFILL.**

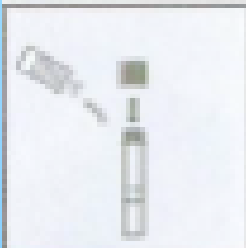


Replace the funnel with the fluid cap.

Remove the funnel from the tube. Screw on the enclosed cap **TIGHTLY** to release the solution that will stabilize the DNA in your saliva.



Tighten to release stabilizing fluid.

You will know it works when the blue solution from the cap has emptied into the tube.

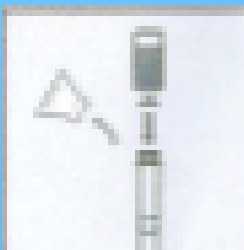

Replace the empty fluid cap with the dead-end cap.

Screw on the dead-end cap **TIGHTLY** to seal your saliva and stabilizing fluid in the tube.



Shake the tube for at least five seconds.

This will ensure your sample mixes thoroughly with the stabilizing solution.

This kit is designed for the collection of human saliva samples.

Cap choking hazard. Keep out of reach of children. Wash with water if liquid solution comes in contact with eyes or skin. **DO NOT INGEST LIQUID.**

Made in USA, Patent Pending
For research use only. Not for use in diagnostic procedures.